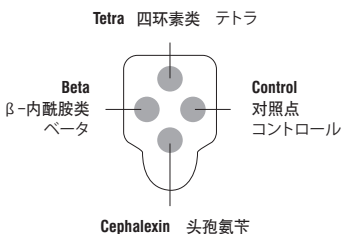
## Results

Resultados • Resultados • Résultats • Wyniki • Ergebnisse • Risultati • 結果 • 結果

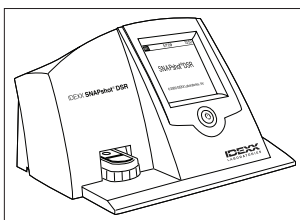
	Beta	Tetra	Cephalexin
	⊖	⊖	⊖
	⊕	⊕	⊕
	⊕	⊖	⊕
	⊖	⊕	⊖
	⊕	⊕	⊖
	⊖	⊖	⊕

OR

O OU ODER 或  
OU LUB OPPURE または

### IDEXX SNAPshot\* DSR Reader

**Beta = sample spot 1**

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Beta = ponto de amostra 1

Beta = spot échantillon 1

Beta = punkt próbki 1

Beta = Probenpunkt 1

Beta = punto del campione 1

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**It is recommended that positive samples be further analyzed. Cephalexin is a member of the beta-lactam class of antibiotics. A positive cephalosporin result should be treated as a positive beta-lactam result.**

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\*SNAP, SNAPduo, SNAPshot and Simplicity from IDEXX are trademarks or registered trademarks of IDEXX Laboratories, Inc. or its affiliates in the United States and/or other countries.

Patent information: [idexx.com/patents](http://idexx.com/patents).

© 2016 IDEXX Laboratories, Inc. All rights reserved. • 06-0001897-01

**IDEXX**

Simplicity from IDEXX\*

## IDEXX SNAPduo\* ST Plus Test

The SNAPduo\* ST Plus is a screening test validated to detect beta-lactam and tetracycline residues in commingled cow milk in accordance with European regulations. Review the sensitivity table on the SNAPduo\* ST Plus product page on [idexx.com/dairy](http://idexx.com/dairy) for detection levels of individual drug residues. If you have any questions, please call IDEXX Technical Services.

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### Technical assistance • [idexx.com/dairy](http://idexx.com/dairy)

US/Canada/Latinoamerica	+1 800 321 0207	日本	+021 6127 9528
EU	00800 727 43399	Within AU	+1300 44 33 99
中国	+010 84423133	Within NZ	+0800 102 084

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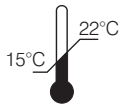
## Ensure test and sample are at correct temperature before use.

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
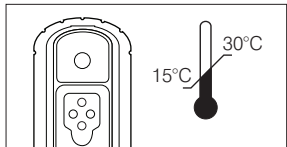
15°C 30°C

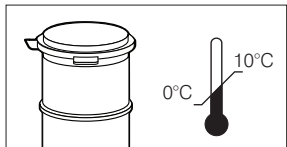0°C 10°C

**Precautions:** Thoroughly mix milk before testing. Do not use abnormal-looking milk or expired tests; do not mix components from different kits. If the control spot fails to develop color, retest the sample.

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### 1 Transfer milk and swirl until dissolved, but no more than 15 seconds.

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